


Hybridoma growth on a 10% XerumFree enriched medium compared with a 10% FBS enriched medium.

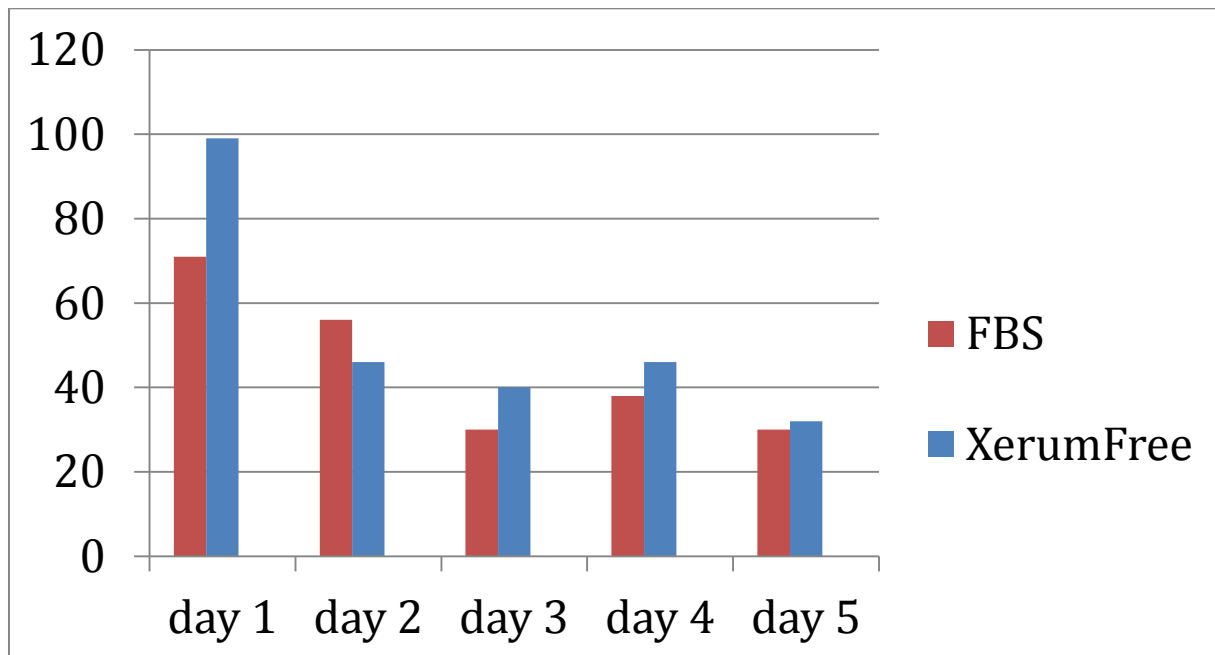

Growth rate of hybridoma cells in %

Every day at PM, cells were passaged in a 5/15 dilution.
This table shows the total cells at the time of the countings

	# of cells/mL (10 ⁶)		% Growth rate FBS	% Growth rate TNC Bio
	10%FBS	10% TNC Bio		
Day 0 PM	0.24	0.01		
Day 1 AM	0.85	0.92	71.37%	98.94%
Day 1 PM	0.37	0.40		
Day 2 AM	0.84	0.74	56.35%	45.95%
Day 2 PM	0.43	0.40		
Day 3 AM	0.62	0.67	30.11%	40.30%
Day 3 PM	0.47	0.40		
Day 4 AM	0.75	0.74	37.78%	45.95%
Day 4 PM	0.43	0.43		
Day 5 AM	0.62	0.64	30.11%	32.29%

Results are from a USA based monoclonal antibody manufacturer.